


VIRTUAL DISSECTION

3D Virtual Dissection


Installation Steps – Windows Vista, Internet Explorer

Step 1 – Close out of any other running programs or browser windows.

Step 2 – Click on the Installation Download link:

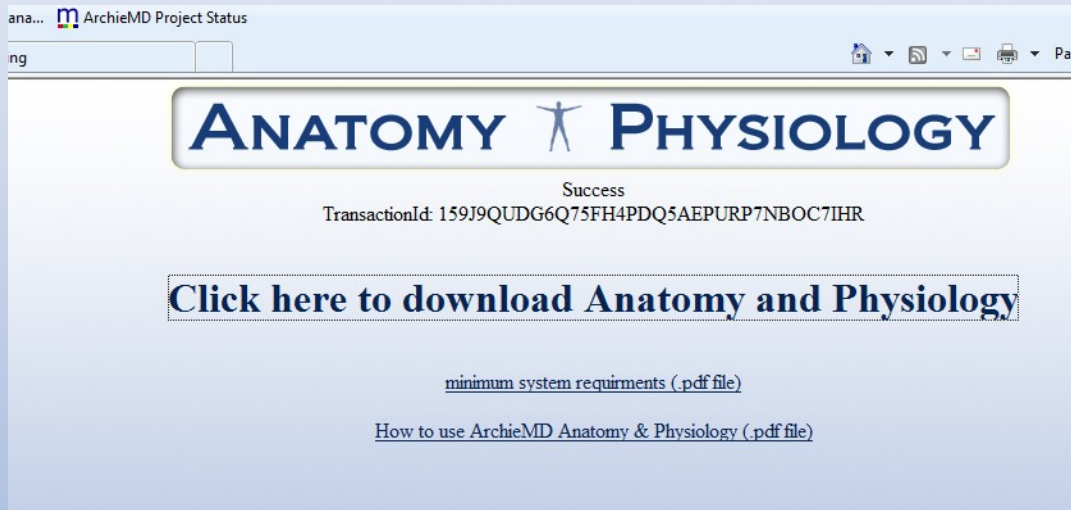

Step 3 – Click on run so that the file will automatically run when the download is complete:

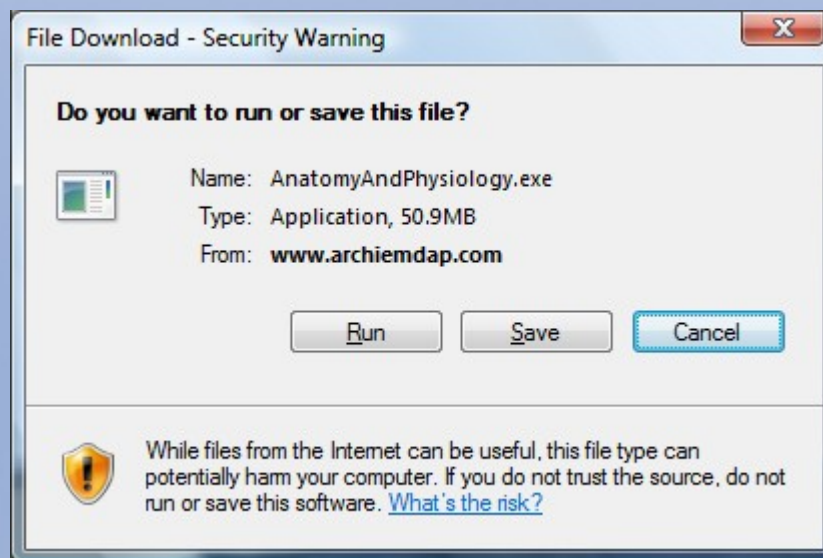


The following warning may appear when the file is finished downloading:




Click "Run".


Once the file finishes downloading, a warning from the User Account Controls may appear. Click "Allow" to proceed with the installation.

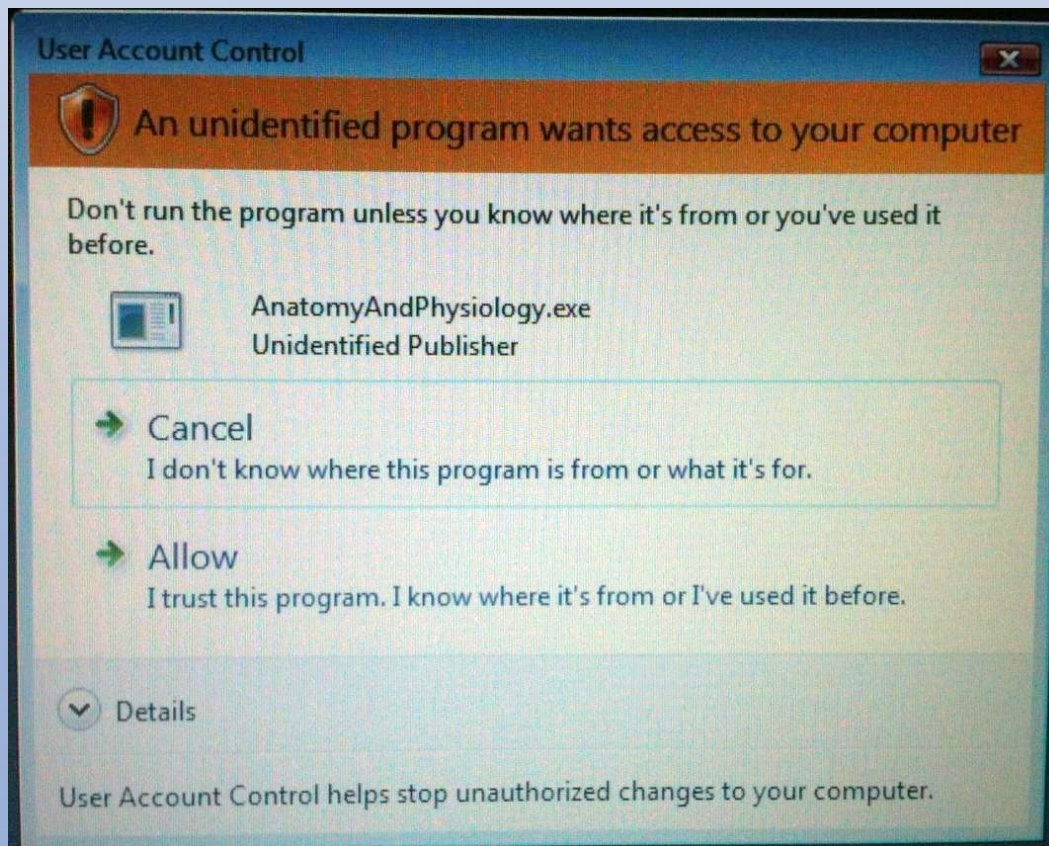


Step 4 – The setup wizard will now launch. Close all other applications and programs before clicking next.

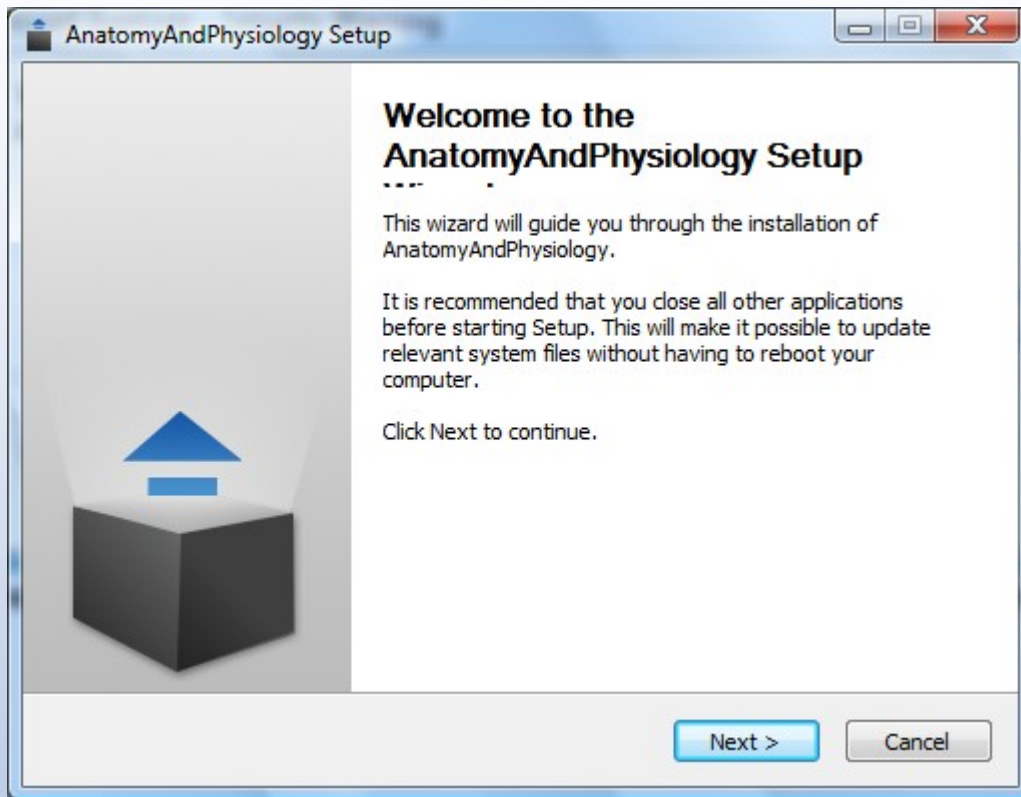

Step 5 – At the next screen, you can decide if you would like the installation wizard to install a desktop shortcut along with Virtual Dissection. This will make the program easier to find once installation is complete. Click “Next” when you are ready to continue.

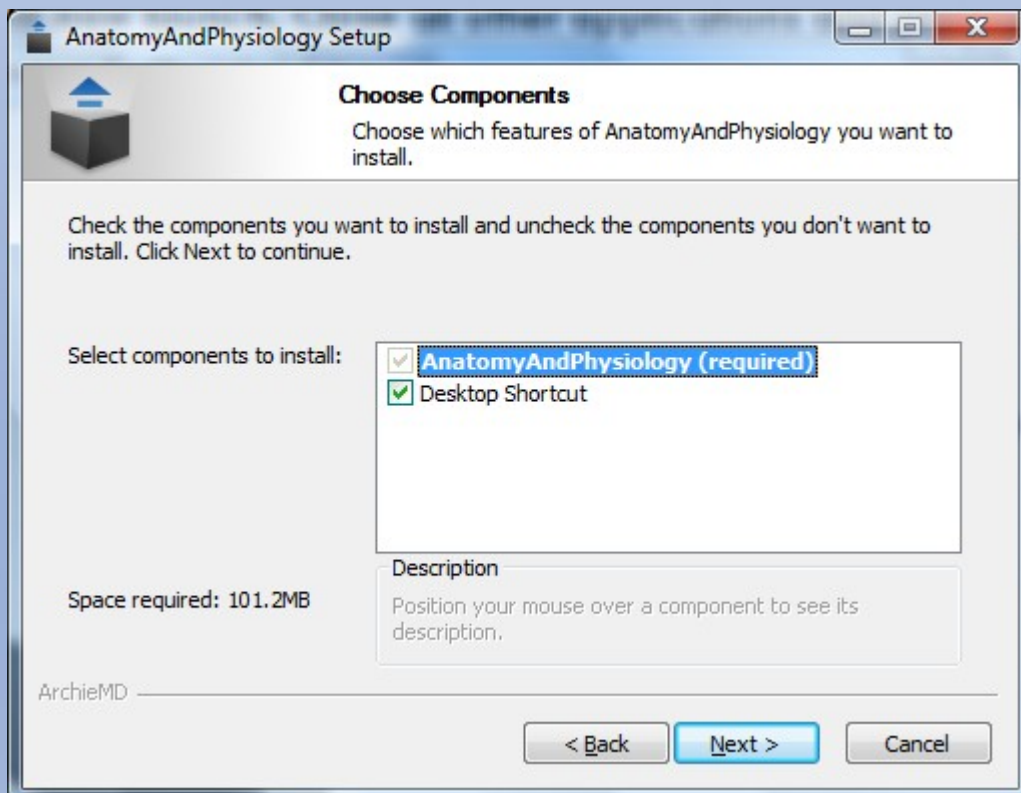


Step 6 – On the next screen, you will choose where Virtual Dissection will install. Choose a location that you will remember and will be able to locate later.

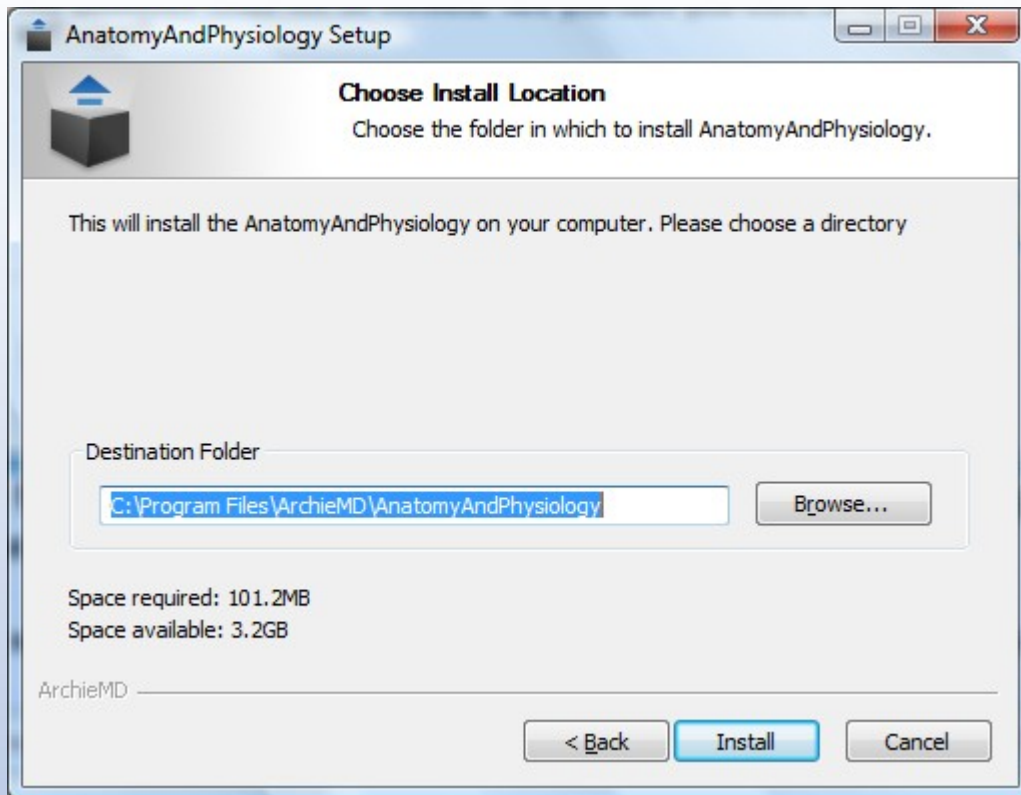


If you would like to change the location Virtual Dissection installs at, simply click “Browse...” A new window will appear, allowing you to choose the appropriate location.

Step 7 – Virtual Dissection will now install. Please wait for it to finish.

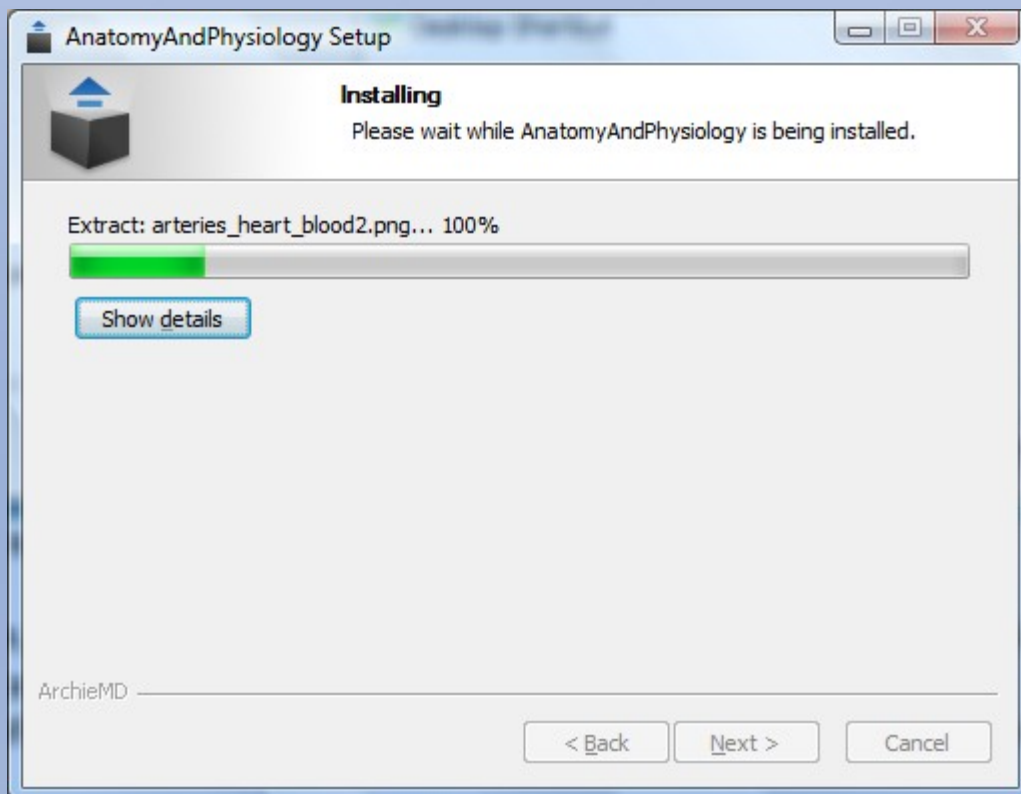

Once the wizard is finished installing Virtual Dissection, you can click “Finish” to close the installer. If you would like to immediately use Virtual Dissection, ensure that the check-box next to “Launch Game Now” is enabled.

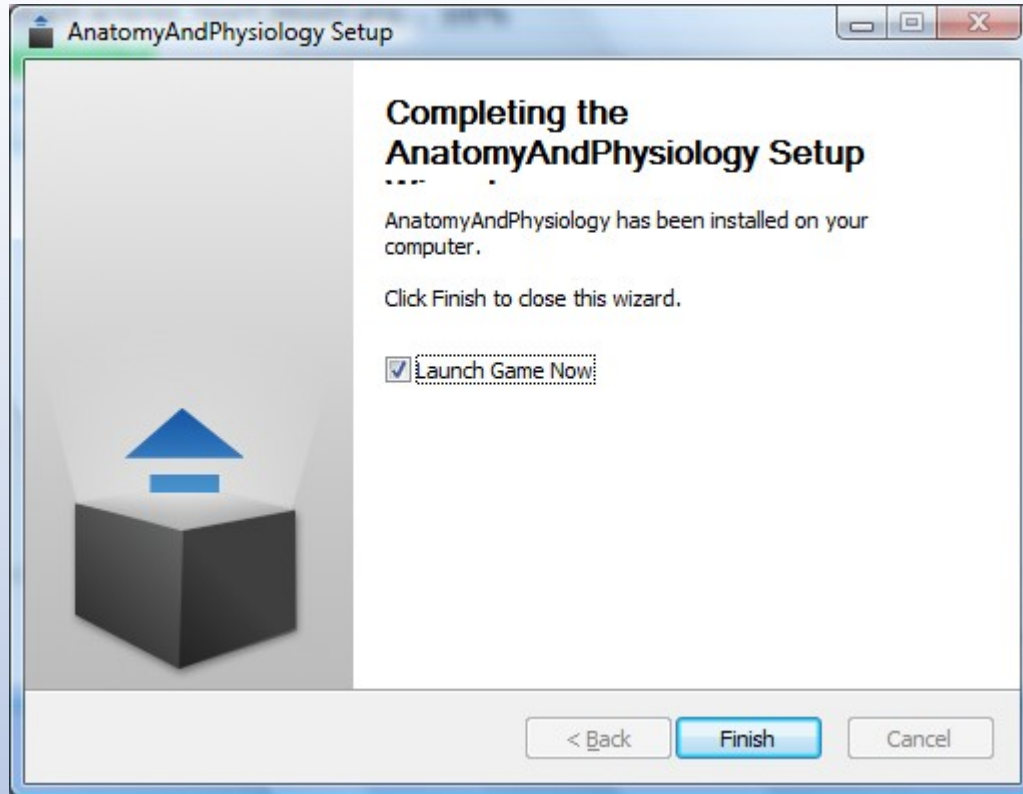

That's all! Enjoy Virtual Dissection!